



MicroSight Focal Tip Instructions

Included with your microscope are five focal tips which allow for easy magnification at specific points. Your microscope will focus anywhere between 10x and 200x and that is a very large range of magnification.

When the specimen is located at the bottom rim of the focal tip, the pre-calibrated tips will magnify at the labeled magnification: 15x, 30x, 50x, 150x, and 100x/200x. (The 5MP MicroSight does not use the 150x focal tip and has only four focal tips.)

Simply pop off the curved lens that has two focal points—100x/200x—and replace it with one of the other tips we provide you and then refocus. Don't be afraid to touch the focal tip down to the surface of your specimen. If your specimen is large or three-dimensional, simply angle the focal tip to one side of your specimen nearest the rim of the tip and then refocus.

When you are using the 100x/200x focal tip, place the focus wheel facing you and turn the focus wheel all the way to the right and then slowly turn the wheel to the left. The low magnification will come into view first at 100x, and when you keep turning past that point, you will soon reach the high magnification point, 200x.

Have fun exploring the hidden world!
Carol

Southern Science Supply
1248 Austin Highway #106
San Antonio, TX 78209

Toll-free: 877-968- 7522
sales@southernsciencesupply.com

www.southernsciencesupply.com